

PROTECTIVE PUMP ENCLOSURE

PART NUMBER #510000

DESIGNED FOR WATSON MARLOW PUMP 120V
(PUMP SOLD SEPARATELY)



**PLEASE FOLLOW THESE GUIDELINES
FOR USE OF THE PROTECTIVE PUMP
ENCLOSURE WITH YOUR WATSON
MARLOW PUMP**



Rotate spring loaded door closure screw to the left. Unscrew lid. Pull lid down.



Cut grommet.



Insert Power Cord through the hole in the back of the enclosure.



Slide black grommet over electrical cord. Grommet needs to have slit in it.



Plug the power cord into the pump.



Slide pump into the case.



Place silicone section of the line set into clamshell, close clamshell over tubing. Push pump securely to the back.



Align tubing with cut out grooves of sidewalls of the enclosure. Close door and verify tube is not being pinched. Secure the door.



**IF USED PROPERLY, THE PUMP ENCLOSURE WILL SIGNIFICANTLY
EXTEND THE LIFE OF YOUR PUMP**

