

# PORTABLE PERISTALTIC PUMP

Part number: 500200



## USAGE GUIDE FOR THE OPERATION OF THE PERISTALTIC PUMP

On/Off Button

Pump Speed Control

Peristaltic Pump

Pump Speed Display

Open clamshell. Place soft section of the tube along rollers. Close clamshell over the tube.

### Tips and Tricks for Ease of Start up and Use:

1. Charge the unit prior to using. Charging will take about 2 hours. A full charge will last up to 11 hours.
2. Press Green On/Off Button to start (Button will illuminate when on).
3. Turn the Pump Speed control dial clockwise (to speed up) and counterclockwise (to slow down).

Speed control turned all the way to 100% equals 90ml per min. 50% equals 45ml per min. Please see sticker on top of the unit for further power setting details.

4. When sampling is complete, press the green button to turn the unit off.

Pump Motor

Battery Pack

On/Off Switch



Battery Charging Cord